

## Commissioning of our Mini-Hydrogen electrode

Before you measure, you must insert the supplied hydrogen cell (button cell) into the electrode head and place the mini hydrogen electrode in water for 24 hours preferably.

**For the functioning of the electrode it is important to stick to the following instructions:**



Unscrew the black cap.



Take the replacement hydrogen cell and remove the protective foil. There are two openings visible. If any adhesive residue remains, remove it by rubbing with a cloth.



Insert the hydrogen cell with these two openings down into the recess.



Turn the cap back in.



Mark month and year on the label. Attach the sticker to the electrode head.



Remove the protective cap from the electrode carefully.



Place Mini-HydroFlex in water for 24 hours.