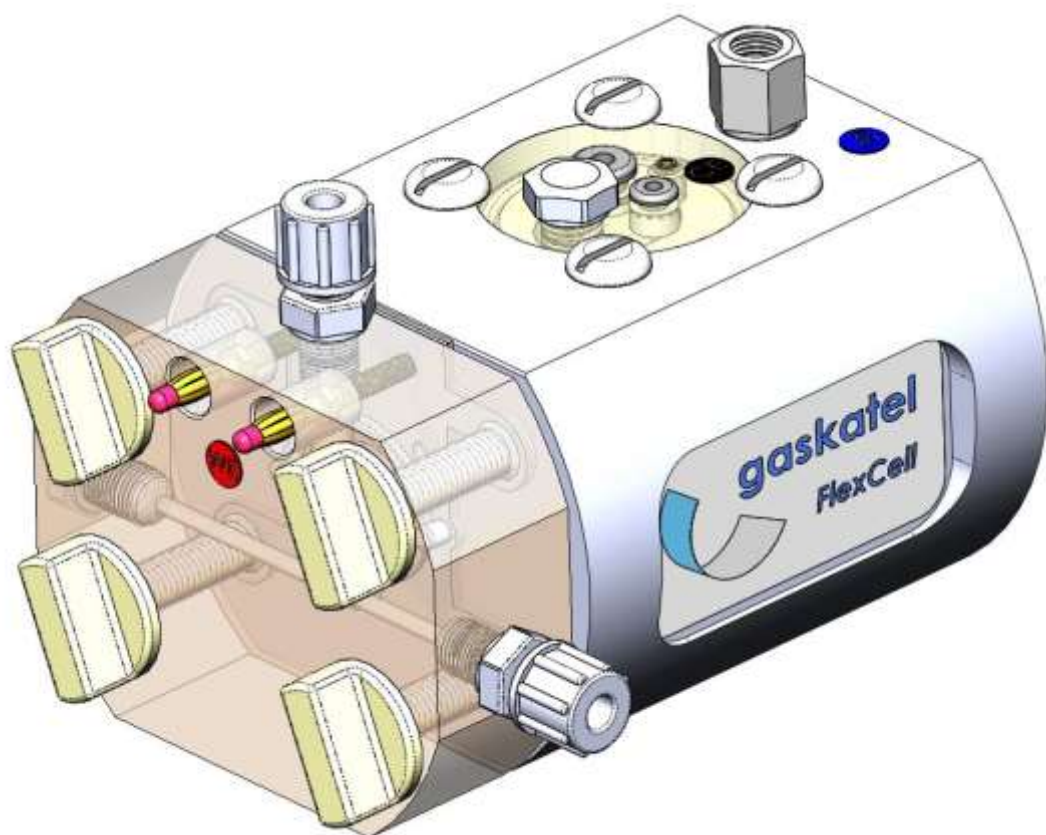
Test Cell FlexCell

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Operating Instructions

Table of contents

Overview	2
Delivery scope	2
Declaration of conformity	4
General.....	4
About these instructions	4
Intended use	4
Safety instructions.....	5
Use	6
Device overview.....	6
Operation	12
Operating conditions	17
Technical data.....	17
Disposal.....	18
Exclusion of liability	19

Overview

Delivery scope FlexCell PTFE 83100

- 1 FlexCell PTFE with
 - 1 Working electrode holder PSU (80547-PSU)
 - 2 Straight connections ECTFE G1/8" (80013)
 - 2 Multilam plugs (89030 (89030)
 - 3 PTFE Pins Ø 4 x 9.5 mm (80029)
 - 1 Screw plug G1/8" PTFE (80031)
 - 1 Counter electrode compartment PTFE
 - 1 PSU-cover (80515)
 - 1 Grommet EPDM 4/8/11-3 (80040)
 - 1 Grommet EPDM 3/5/8-2 (80041)
 - 4 Round head screw slotted M6x16 PA 6.6 (83004)
 - 1 O-ring EPDM, 40 x 3 mm (80541)
 - 1 Screw plug G1/8" PTFE (80031)
 - 1 Spiral counter electrode Pt-Ir, Ø 0.25 mm (830088-v02)
 - 3 Nafion tubes (83010)
 - 2 Silicone sealings (83001)
 - 4 Wing screws M8x50 PA6.6 (83006)
 - 2 Heating elements 200W, 30V cast in silicone rubber (83013)
 - 1 Screw plug G1/4" PP (80030)
 - 1 Screw plug G1/8" PTFE (80031)
 - 1 Connection cable for counter electrode CE (89013)
 - 2 Connection cables for working electrode WE (89017)
 - 4 Connection coupling red (89022)
 - 1 Connection coupling black (89023)
 - 2 O-Rings EPDM, 24 x 2 mm (80321)
 - 1 Reducing nipple G1/8" G1/4" PP (80017)
 - 1 Socket wrench screwdriver 6 x125mm 6kt.
 - 1 Manual

Delivery scope FlexCell PP 83200

- 1 FlexCell PP with
 - 1 Working electrode holder PMMA (80547-PMMA)
 - 2 Straight connections PP G1/8" (80012)
 - 2 Multilam plugs (89030 (89030)
 - 3 PTFE Pins Ø 4 x 9.5 mm (80029)
 - 1 Screw plug G1/8" PP (80003)
 - 1 Counter electrode compartment PTFE
 - 1 PSU-cover (80515)
 - 1 Grommet EPDM 4/8/11-3 (80040)
 - 1 Grommet EPDM 3/5/8-2 (80041)
 - 4 Round head screw slotted M6x16 PA 6.6 (83004)
 - 1 O-ring EPDM, 40 x 3 mm (80541)
 - 1 Screw plug G1/8" PP (80003)
 - 1 Spiral counter electrode Pt-Ir, Ø 0.25 mm (830088-v02)
 - 3 Nafion tubes (83010)
 - 2 Silicone sealings (83001)
 - 4 Wing screws M8x50 PA6.6 (83006)
 - 2 Heating elements 80W, 30V cast in silicone rubber (83012)
 - 1 Screw plug G1/4" PP (80030)
 - 1 Screw plug G1/8" PP (80003)
 - 1 Connection cable for counter electrode CE (89013)
 - 2 Connection cables for working electrode WE (89017)
 - 4 Connection coupling red (89022)
 - 1 Connection coupling black (89023)
 - 2 O-Rings EPDM, 24 x 2 mm (80321)
 - 1 Reducing nipple G1/8" G1/4" PP (80017)
 - 1 Socket wrench screwdriver 6 x125mm 6kt.
 - 1 Manual

Optionally available accessories

Hydrogen reference electrode Mini-HydroFlex 81020

Temperature control box 83900

Analyte compartment PTFE 83110

Analyte compartment PP 83210

Declaration of conformity

 This product complies with all EU directives applicable to this product.

General

About these instructions

Read these instructions carefully. To ensure a long service life and reliable use, all instructions mentioned in this manual must be observed and followed. Please keep these instructions handy.

You can access the instructions and further information at www.gaskatel.com.

Intended use

The FlexCell test cell is suitable for electrochemical measurements on gas diffusion electrodes and other flat samples such as metals.

We supply appropriately cut silicone seals and O-rings made from EPDM. Make sure that the seals are well seated before assembling the test cells.

The test cells are designed for the typical three-electrode set-up. A counter electrode is already integrated.

Both test cells can be heated via integrated heating elements. They may only be operated with the [Gaskatel temperature control box 83900](#).

They are connected in parallel to the temperature control box.

The [test cell FlexCell PTFE](#) may be operated up to **150°C**.

The test cell [FlexCell PP](#) measuring cell may be operated up to **80°C**.

The [filling level of the electrolyte must be checked regularly](#), as the level can change significantly during the measurement.

The measuring cells may only be operated in a laboratory environment.

Safety instructions

The FlexCell test cells are only intended for the above-mentioned purpose.

The manufacturer is not liable for damage resulting from improper use.

Observe the following instructions, otherwise the test cells may be damaged or measurement results may be falsified.

When working with chemicals, all relevant safety regulations of the manufacturer and the laboratory must be observed.

The cell heater and / or the cell current you apply can heat the test cell over 80°C/150°C.

The test cells must be placed on a base that is insensitive to heat. Do not place heat sensitive things on the half cell.

Do not place inflammable things near the half cell.

Pay attention to a correct lining of your measurement and heating cable.

Select cable insulations with a reasonable temperature specification. Else there is a risk of an electrical shortcut and fire hazard.

Electrolyte vapours are constantly emitted from the test cell according to the vapour pressure curve.

The test cell should be operated on a laboratory tray that can safely hold the entire filled electrolyte volume in case of leakage.

Protect electrical / electronic equipment by a large distance, partitions or operation under a fume cupboard.

Wear safety glasses.

The electrolyte can splash out of the test cell in the event of pressure surges on the gas inlet or when the boiling point is reached.

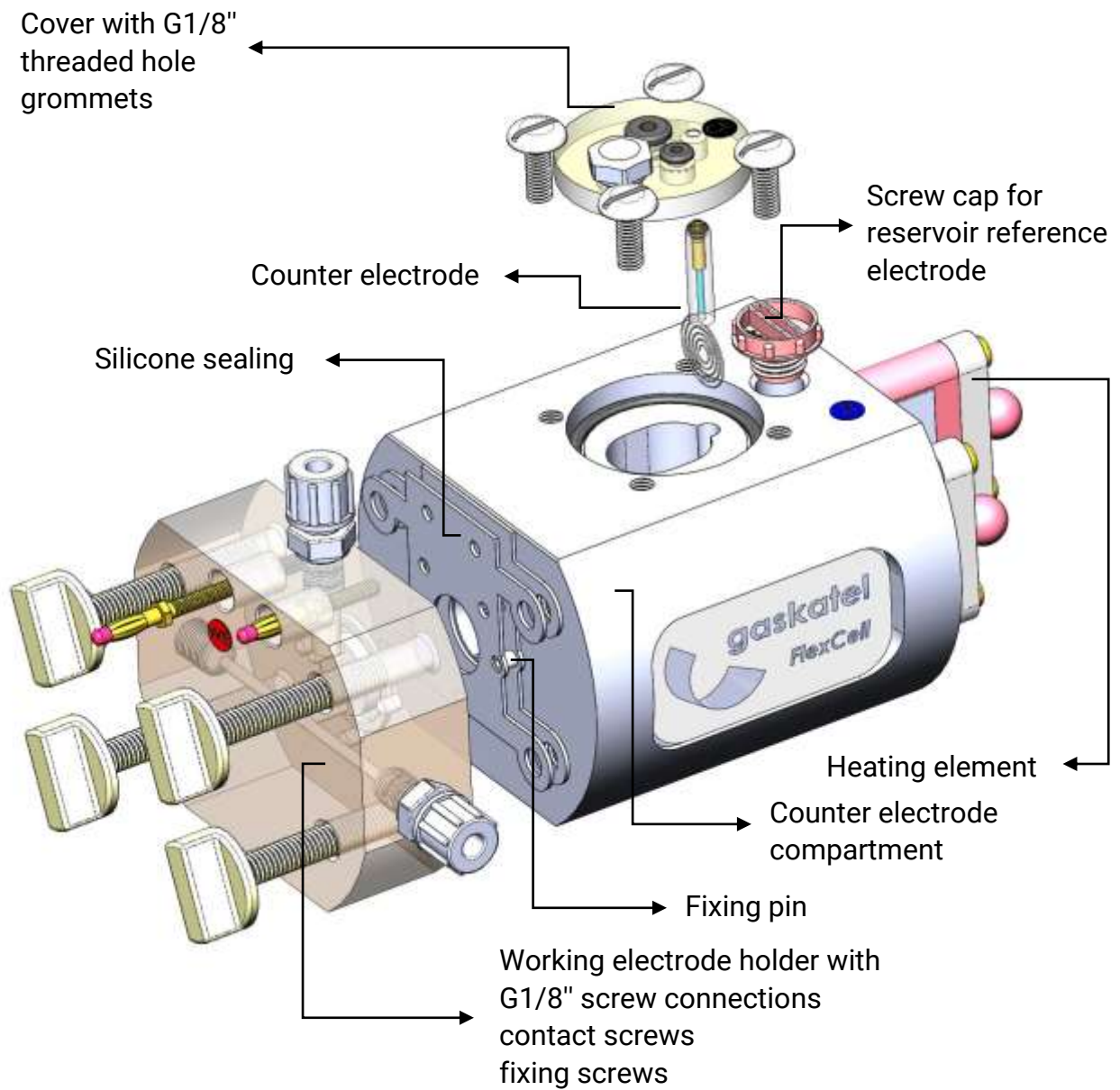
The operator of the test cell must be equipped with adequate laboratory protective equipment according to the safety data sheets and specifications for your electrolyte.

Use


Device overview

Layout

Principle of test cell FlexCell

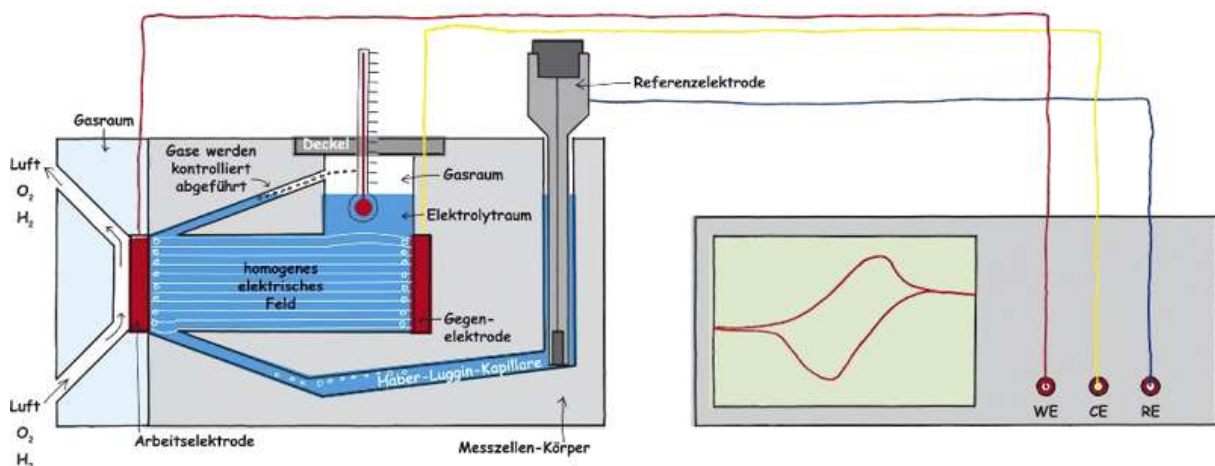To characterise electrochemical processes, the so-called three-electrode arrangement is used. The three-electrode arrangement consists of the working electrode (measuring electrode), the reference electrode (sensor electrode) and the counter electrode. These kinds of measurements are usually controlled by potentiostats or galvanostats. Electrochemical test cells have not been standardized yet. Therefore, there are many different types, some with serious errors. Through many years of experience, especially in the fuel cell and battery sector as well as corrosion measurements, Gaskatel has successfully constructed an electrochemical test cell that largely eliminates or minimises these known errors.

Made of plastic

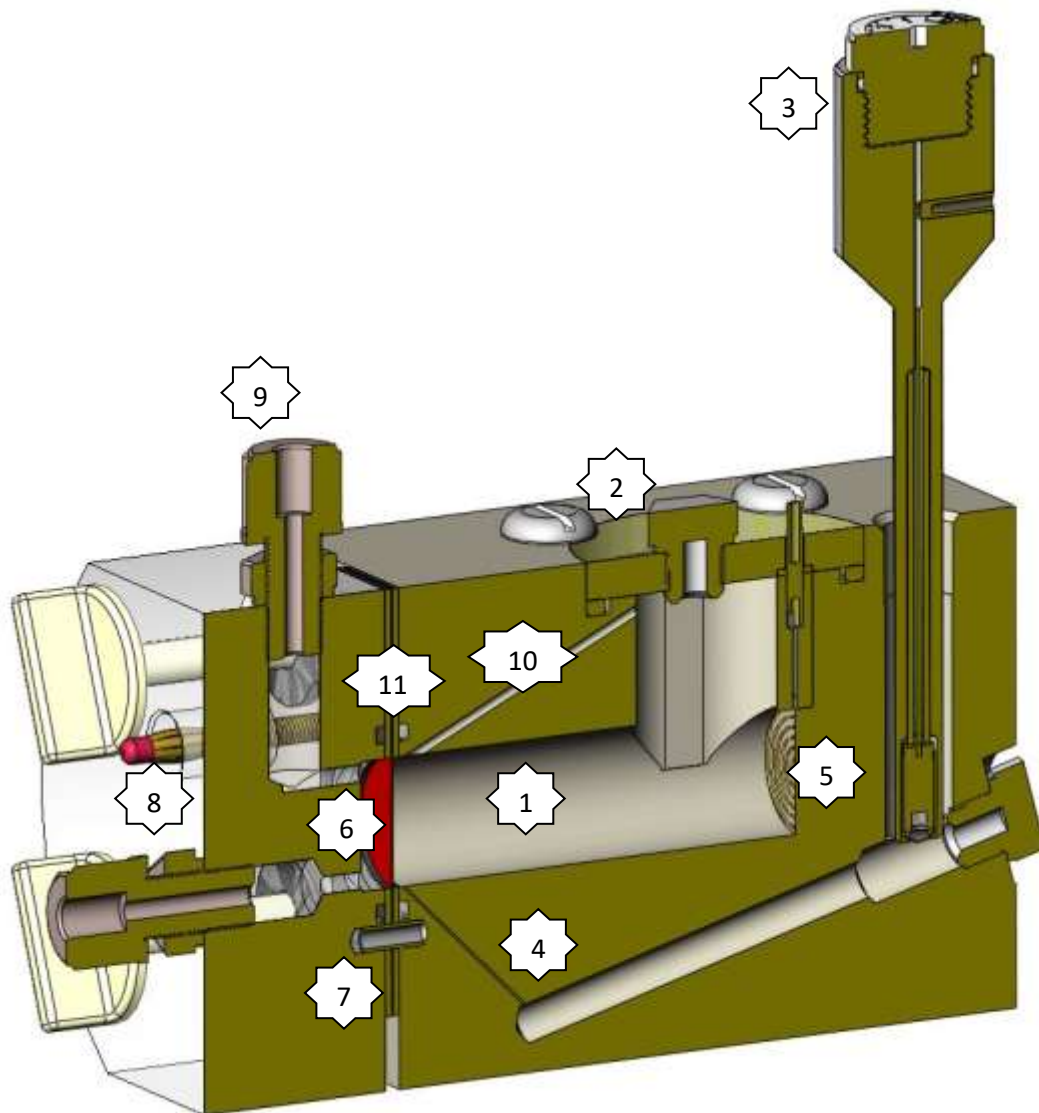
In contrast to the usual measuring cells, Gaskatel test cells are made of PTFE or PP, because glass is a poor material. For one thing, it breaks quickly. For another, it is chemically stable only to a limited extent, because silicates dissolve. This is a particular problem with aggressive solutions such as alkalis.

But also, for light sensitive electrolytes glass is an insufficient material.

In very poorly conductive electrolytes, the conductive silicate layers cause an uneven field line distribution.

Gaskatel test cells made of PTFE and PP are characterised by high mechanical stability and high chemical resistance. The plastics have no influence on the field line distribution.

Detailed view of the test cell FlexCell

Counter electrode compartment (1)

Only test cells with a tubular arrangement with equally sized measuring and counter electrodes guarantee a parallel field line distribution.

Parallel field lines in turn guarantee parallel equipotential surfaces and thus a reproducible potential measurement, as the reference electrode measures on these potential surfaces.

Gaskatel test cells have such parallel field lines, as the working and counter electrodes have the same size and are arranged parallel to each other due to the tubular electrolyte space.

To minimise water vapour losses, the cell is closed with a cover made of PSU (2).

Reference electrode (3)

Gaskatel test cells are optimised for the use of the reversible hydrogen reference electrodes (RHE) Mini-Hydroflex.

In contrast to the usually used calomel or silver electrodes, these hydrogen reference electrodes do not introduce any foreign electrolyte into the measuring solution. This is very important for corrosion measurements, for example, where the smallest amounts of chloride can drastically influence and falsify the results.

Most electrochemical reactions are pH dependent. This is often compensated by the use of Mini-HydroFlex, because Mini-HydroFlex itself is also pH-dependent.

Position of the reference electrode – Haber-Luggin-capillary (4)

For precise potential measurements it is necessary that the reference electrode is positioned close to the working electrode.

To reduce the voltage drop across the electrolyte, the so-called IR drop, Haber-Luggin capillaries are used. However, for measurement results to be comparable between cells, the Haber-Luggin capillary must meet very precise and reproducible requirements.

The field lines in the electrolyte must not be disturbed (small diameter). In turn, a large diameter is necessary for low electrolyte resistance. Therefore, the capillaries in Gaskatel test cells are located outside the electrolyte space. This allows a slightly larger diameter without disturbing the field lines.

The capillary must end at a defined, fixed distance from the working electrode. From a quality assurance point of view, it is also important that all these factors are kept constant for different half-cells.

This is where the precision of CNC manufacturing in plastic helps. The Gaskatel measuring cells have very precise Haber-Luggin capillaries that end at the working electrode. This makes the measurements very precise and comparable.

The capillaries must not be blocked by gas bubbles, as this leads to contact losses during the potential measurement. The resistance then increases dramatically and the potentiostat starts to oscillate.

Therefore, the Haber-Luggin capillaries in the FlexCell measuring cells are filled with a solid electrolyte.

Counter electrode (5)

The counter electrode must be electrochemically resistant so that no foreign ions are introduced into the electrolyte.

The overvoltage should be low, otherwise the temperature in the electrolyte will rise sharply.

The FlexCell test cells have a platinum-iridium spiral with a wire thickness of 0.25 mm as a counter electrode. The wire is connected to a gold-plated 2 mm banana socket on the cover.

Working electrode (6)

The working electrode is placed between two seals. To prevent the seals from slipping, there are 3 PTFE pins (7). For fitting between these pins, the sample must be cut to the dimensions 30 mm x 50 mm.

If you want to use electrodes of other geometries, you can simply remove these PTFE pins.

If the working electrode is a gas diffusion electrode with a non-conductive gas diffusion layer such as porous PTFE, then the PTFE film must be removed about 1 cm from the edge so that the sample can be contacted there. The catalyst mass can also be removed carefully on this piece. This prevents air from entering the sample from the outside and is especially important for measurements on nickel hydrogen electrodes.

Contact working electrode (8)

Two 4 mm gold-plated multilam plugs press onto the working electrode from the rear. If the contact is not sufficient, the multilam plugs can be tightened with a socket spanner. Make sure that these plugs are clean. Replace these plugs if the gold-plated layer no longer looks sufficiently good. This also applies to all measuring cables.

Gas supply (9)

For measurements on gas diffusion electrodes, these must be supplied with gas. Here, it is done via the working electrode holder, made of transparent plastic so that the electrode can be observed. It has two connections: the gas inlet at the top and the gas outlet at the bottom. If liquid gets through the electrodes, it can be removed with the gas flow.

Gas bubbles (10)

There cannot be any electrolyte where the gas bubbles stick to the electrode. Gas bubbles can block the reaction surface. Local elements can form. Gaskatel test cells have an additional opening through which any gas bubbles that form are removed in a controlled manner.

Crevice corrosion (11)

If the sample is not properly sealed, an electrochemical reaction may take place underneath the sealing or even outside of the active area. A proper sealing is essential. Oxygen may only enter in a controlled manner.

Flat sealings have proven successful with very noble materials, thin samples or gas diffusion electrodes.

O-ring seals should be used with base materials such as steel, aluminium, etc.

The measuring cells are therefore supplied with the appropriate O-rings and appropriately cut silicone flat sealings.

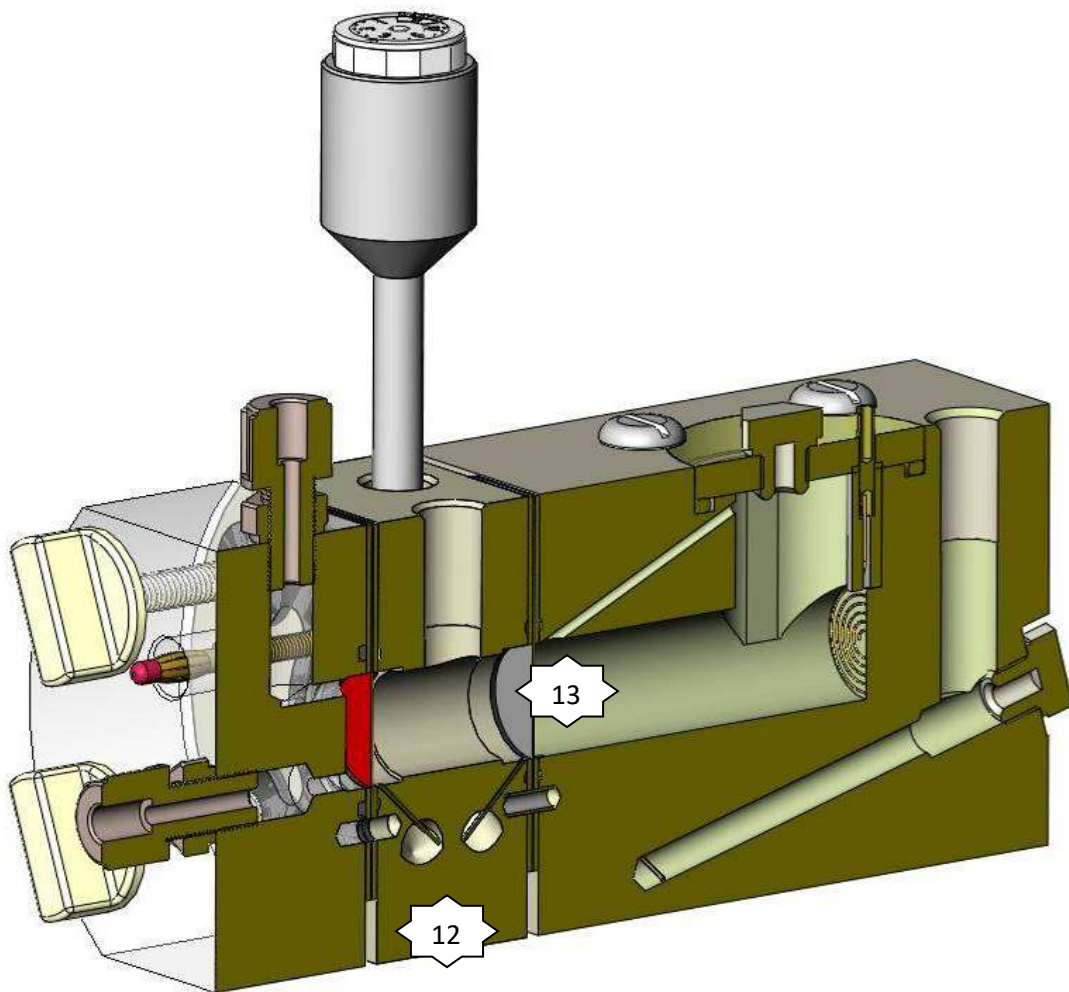
Reaction products of the counter electrode – analyte compartment (12)

Often very aggressive ions are produced at the counter electrode. Depending on the electrolyte, peroxides, perchlorates or persulphates etc. can be produced. Even in the smallest quantities may initiate corrosion on the working electrode.

Therefore, Gaskatel offers an intermediate plate (analyte compartment) to place a membrane (13) between the counter and working electrode. In this way, working and counter electrode are separated from each other.

The reference electrode must then be installed in this analyte compartment.


Corresponding Haber-Luggin capillaries with solid electrolyte are of course available in the analyte compartment.

Temperature

Many processes take place at higher temperatures. Therefore, it is often necessary to heat the test cells.

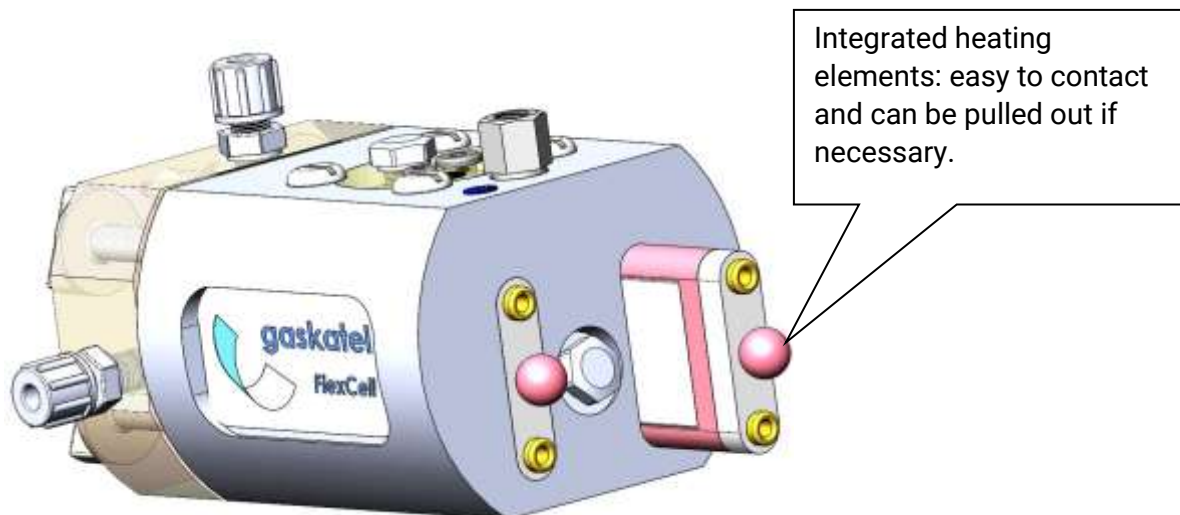Gaskatel FlexCell test cells can be heated via integrated heating elements. These are placed into corresponding boreholes and can simply be pulled out before cleaning the test cells.

Please use our temperature control box 83900 for temperature control, because only a well-regulated temperature ensures comparable measurement results.

Operation

1. Assembling the test cell

First, the working electrode may have to be cut to size:

Width 30 mm, length 50 mm

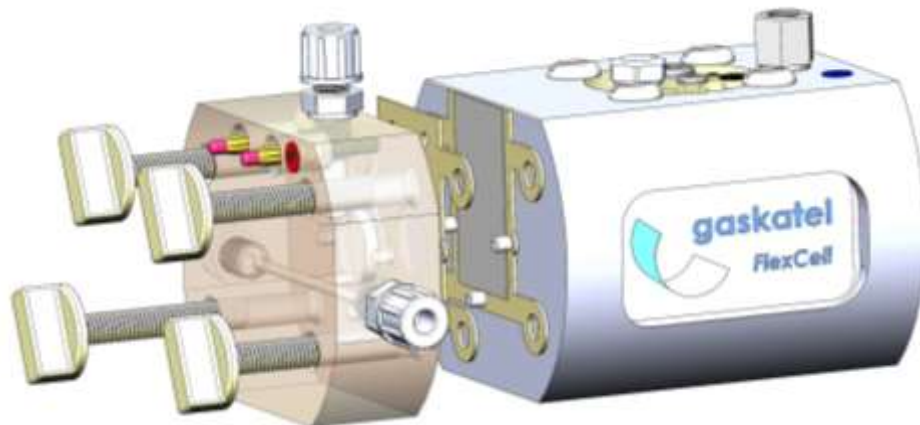
The thickness of the working electrode is almost arbitrary.

The active area in the electrolyte is 3 cm².


The contact screws in the working electrode holder must be turned out a little so that the electrode cannot touch these contacts at first.

The cut working electrode is placed between the two silicone sealings or EPDM O-rings.

The fixing pins ensure that the electrode cannot slip.

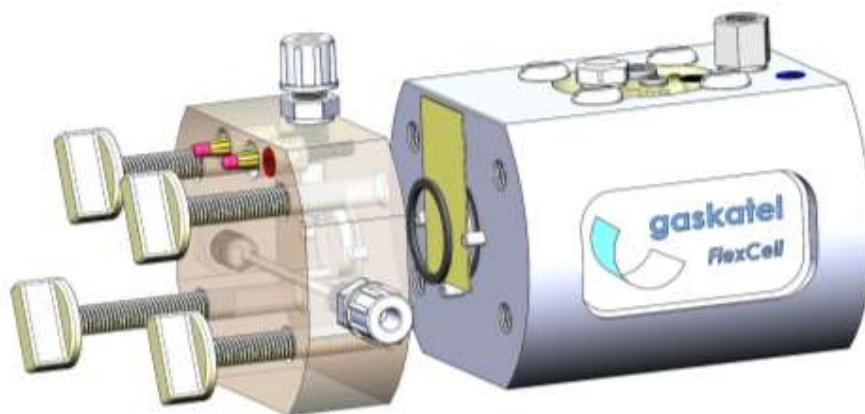FlexCell with inserted gas diffusion electrode between the silicone sealings

FlexCell with inserted metal sheet between the O-rings

When screwing the cell components together, make sure that the screws are tightened diagonally to each other one after the other. There must be no gap between the components. Tighten the screws only hand-tight, especially when using the flat gaskets. Leakage is usually caused by unevenly tightened screws and not by too weakly tightened ones.

If this is the case, you can now screw the contact screws back in with the socket spanner and hand-tighten them so that you contact the electrode.

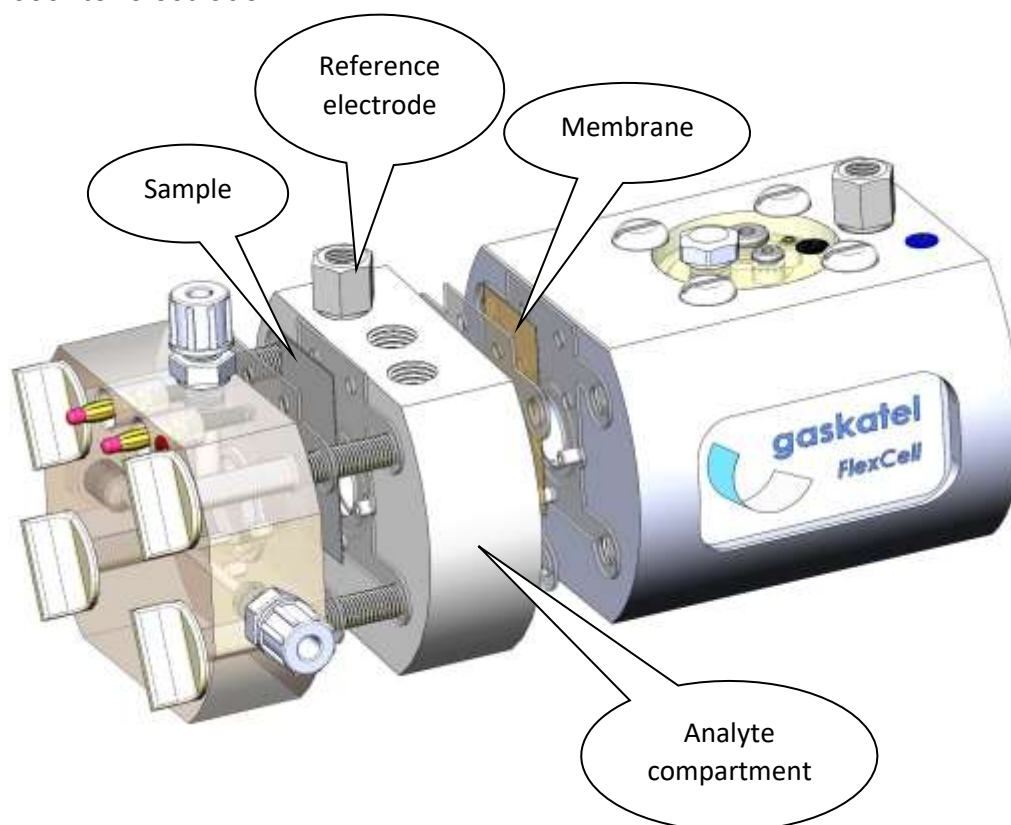
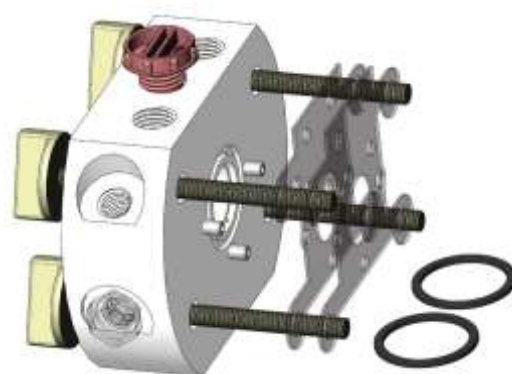
2. Filling of the test cell

The cell can be filled with approx. 30 ml of electrolyte through the opening in the cover. Pour in the last ml slowly; observe the filling level in the hole for the reference electrode, which may rise with a delay. This should be filled to a maximum of halfway to avoid overflowing when setting the reference electrode.

If the hole does not fill with electrolyte, an air bubble in the Haber-Luggin capillary blocks the electrolyte flow. Set the measuring cell at a slight angle so that the air bubbles can escape.

3. Use of the analyte compartment with membrane

In addition, an intermediate plate (analyte compartment) can be inserted. These are optionally available in PP or PTFE and are supplied with the necessary longer screws. A membrane is inserted between the analyte compartment and the counter electrolyte compartment. This membrane separates anolyte and catholyte. The working electrode is thus protected from the reaction products of the counter electrode.

When selecting the membrane, pay attention to the electrolyte resistance and the ion-exchanging properties of the membrane.

For the potential measurement of the working electrode, the reference electrode must be placed into the analyte compartment.

If you have a bad conducting electrolyte, the potentiostat may start to oscillate due to these conductivity problems in the Haber-Luggin capillary. In this case, if you use the analyte compartment with a much lower resistance Haber-Luggin capillary, the problems should be solved.

You can also measure membrane resistances with this arrangement. Two reference electrodes are then needed. One is inserted into the analyte compartment and the second into the bore of the electrolyte compartment as usual.

4. Cleaning of the test cells

Usually, it is sufficient to rinse the cell thoroughly with water after disassembly.

Deposits can be removed mechanically with a suitable laboratory brush.

Stubborn metal deposits such as silver or copper can be removed with nitric acid.

5. Troubleshooting

Error	Possible causes	Check
Cell temperature does not exceed 80°C (PP) 150°C (PTFE)	Bad contact	Check the connecting leads of the heating elements
	PTC heating element in silicone is defect	Measure the resistance of cold the heating elements: PTFE-Test cell: 5 Ohm PP-Test cell: 3 Ohm
Potentiostat is going in overpotetnial	Bad contacts	Check the connecting leads of the potentiostat
	Pt counter electrode	Remove the counter electrode and check its condition
False potential	Gas bubbles in front of the reference electrode	Move the reference electrode a little up and down
	Gas bubbles in Haber-Luggin capillary	Suck air out of the reference electrode hole with a pipette ball
Noisy signal	Impedance of the reference electrode or Haber-Luggin capillary is too high	Adjust the current range of the potentiostat
		Use the analyte compartment: reference electrode must be set in the analyte compartment
Electrolyte leakage	Plastic wing screws not tightened	Tighten the wing screws
	O-ring or silicone sealing is not placed properly	Check the position of the contact screws and fixing pins.

Operating conditions

	FlexCell PTFE	FlexCell PP
Nominal voltage working- / counter electrode:	< 50 VAC < 75 VDC	< 50 VAC < 75 VDC
Nominal current working- / counter electrode:	max. 3 A	max. 3 A
pH-range:	pH -2 bis pH 16	pH -2 bis pH 16
Temperature:	10°C bis 150°C	10°C bis 80°C
Pressure:	max. 200 mbar	max. 200 mbar
Maximum recommended measurement duration:	24 h	24 h

Technical data

	FlexCell PTFE	FlexCell PP
Dimensions: without analyte compartment, protruding screw connections, screws	132 mm x 100 mm x 75 mm	
Electrolyte volume:	30 ml	
Material of counter electrode compartment:	Polytetrafluorethylene	Polypropylene
Material of cover counter electrode compartment:	Polysulfone	
Material of working electrode holder	Polysulfone	Polymethyl- methacrylat
Active sample area:	3 cm ²	
Working electrode sample size:	30 mm x 5 mm	
Contacts working electrode:	2 multilam plugs 4 mm, gold-plated	
Sealings working electrode:	Silicone sealings, O-Rings EPDM	
Screw fitting gas supply:	ECTFE G 1/8" für 6 / 4 mm tube	PP G1/8" für 6 / 4 mm tube
Heating elements (cast in Silicone):	200W, 30V	80W, 30V
Counter electrode:	Spirale Pt/Ir-Draht, Ø 0,25 mm	
Contact counter electrode:	2 mm banana socket, gold-plated	
Fixing screws:	Polyamide PA 6.6	

Disposal

The test cell must not be disposed of with household waste. Please dispose of the test cell in accordance with local regulations in a separate collection for electrical and electronic equipment.

If sent to us free of charge (sufficiently stamped package), we will dispose of the used test cell from our product range free of charge.

Exclusion of liability

The product is subject to continuous further development. For this reason, it is possible that parts of the instructions, technical data or images in these instructions may differ slightly from the product in front of you. The information in these instructions is only intended to clarify how the product is to be handled.

If you have any questions, please do not hesitate to contact us.

Legal claims based on these instructions cannot be asserted.