

## Replacement of the hydrogen cell in Mini-HydroFlex

With a few simple steps, you can replace the hydrogen cell in Mini-HydroFlex hydrogen electrode.

**For the functioning of the electrode it is important to stick to the following instructions:**



Unscrew the black cap and remove the hydrogen cell.



Grease  
O-ring

In the recess in the electrode head you see a small o-ring. Please grease this, e.g. with silicone grease.



Take the replacement hydrogen cell and remove the protective foil. There are two openings visible. If any adhesive residue remains, remove it by rubbing with a cloth.



Insert the hydrogen cell with these two openings down into the recess.



Turn the cap back in.



Mark month and year on the label. Attach the sticker to the electrode head.



24 h  
Aqua  
dest.

Place Mini-HydroFlex in water for 24 hours.